


ISCUSflex QC Charts Installation Instructions


Download and save the installation file (ISCUSflex QC Charts.zip).



Click on Save

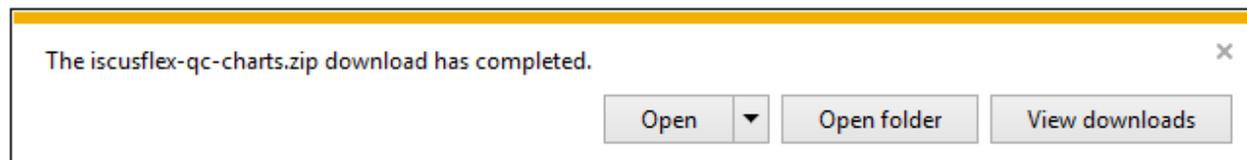

Click on Open Folder

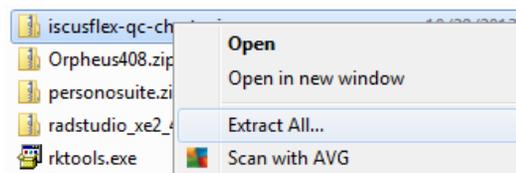

Select Extract All ...

When finished double click on the Setup icon to start the installation.

Name	Date modified	Type	Size
Application Files	10/29/2013 10:32 ...	File folder	
ISCUSflex QC Charts.application	10/29/2013 10:32 ...	ClickOnce Applica...	3 KB
setup.exe	10/29/2013 10:32 ...	Application	425 KB

If .NET 4.5 is not installed it will be automatically installed on the computer. When the installation has finished the software is automatically launched.

An icon is created in the Start menu.

N.B. At first time start of the software you need to enter the license code. The licensed code can be purchased from M Dialysis. A license code is valid for one computer only.